

CANTY

PROCESS TECHNOLOGY

INDUSTRIAL SOLIDSIZER™ Particle Analysis System




ONLINE PARTICLE SIZING


The Canty Industrial SolidSizer™ is an on-line particle sizing tool for lab environments to determine particle size, shape and distribution, thereby eliminating the need for sieve analysis.

THE CANTY ADVANTAGE

The JM Canty Industrial SolidSizer™ is a vision-based sensor used with the Vector System for dry particle size measurement in a production environment as an on-line production tool. The SolidSizer™ includes a B & W camera with a shutter speed control and a variable magnification lens, a light source to generate a uniform light field to display silhouette images of opaque particles, and a vibratory feeder to present sample material to the camera / lights in free fall. The weatherproof camera is equipped with a spray ring that can be charged intermittently to control dust buildup on the outside of the camera view port. The camera video is provided to the Canty Core unit for analysis that extracts particle size information. This sheet identifies the features of the SolidSizer™ and provides ordering information. Typical installations also include the Canty Core unit.

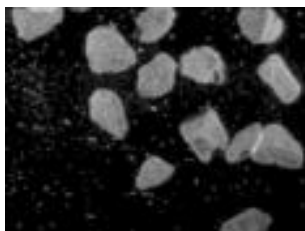

Live Image

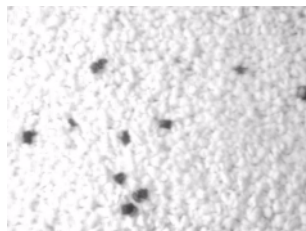

Black Speck Detection

FEATURES

- Particle distribution by major/minor diameter, area, perimeter, etc...
- Real time, on-line particle size analysis
- Adjustable lens magnification (sub micron through boulders – no upper size limit). Consult factory
- ASTM D8200 sieve correlation
- Shape Analysis by length, width, aspect ratio, circularity, solidity etc...
- External video connectors for signal processing by Canty Core unit (optional)
- Optional light filters available for wavelength sensitive materials
- Excel data format available with Canty Core unit

APPLICATIONS

- Pharmaceutical powders
- Polymer pellets and beads
- Agricultural products
- Aggregates
- Crushed rock
- Glass bead applications
- Screen break detection
- Agglomeration control
- Spray dryers and fluid bed applications
- Replaces existing lab screen/sieve systems
- End product Q.C. and lot certification

BENEFITS

- Reduces Lab Screening Labor by 99%
- Output Matches Sieve/Screen Data - Direct Replacement
- Three Month Pay Back

SPECIFICATIONS

- Power: 120V AC / 60 Hz (230V AC / 50 Hz)
- Shutter Speed: variable up to 1/100,000 sec
- Instrument Air: 20 to 30 PSI when required

CANTY
www.jmcanty.com

JM Canty Inc
JM Canty Intl Ltd

Buffalo, NY USA
Dublin, Ireland

Ph: 1 (716) 625 4227
Ph: + 353 (01) 882 9621

Ordering Information

HOW TO ORDER: Select the appropriate symbols to build a part number as shown:

EXAMPLE:

S C - G 2 K 1 2 2 1

VIDEO SYSTEM FORMAT

G - Gigabit Ethernet, B&W, 1600 x 1200

H - Color, 640 x 480

ENVIRONMENTAL RATING

2 - Weatherproof

3 - Explosion Proof

ACCESSORIES

1 - No Accessories

2 - Black Speck Only

3 - Particle Size and Black Speck

INPUT POWER




1 - 120 V AC 60Hz

2 - 230 V AC 50Hz

SYSTEM ILLUMINATION

K - Backlighting

L - Back / Front (Ringlight), Black Speck Only

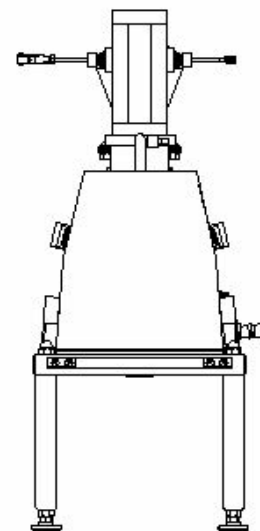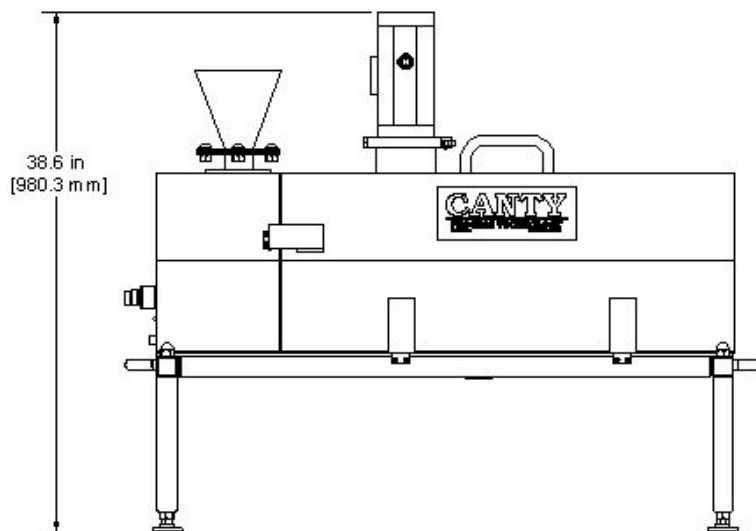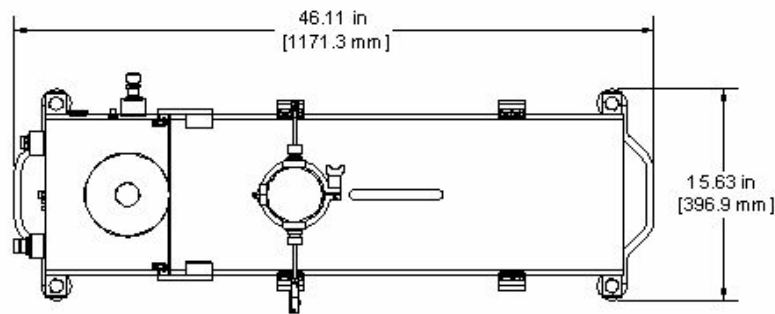

* Fines separator is used to separate small particles (under 80 micron) for optimal particle orientation.
Consult factory if required for your application.

CANTY

www.jmcanty.com

JM Canty Inc
JM Canty Intl Ltd

Buffalo, NY USA
Dublin, Ireland

Ph: 1 (716) 625 4227
Ph: + 353 (01) 882 9621